



## Water determination in dimethyl sulfoxide (DMSO)

HYDRANAL™ Laboratory Report L 141

It is not possible to reproduce water determination in DMSO because it alters the stoichiometry of the Karl Fischer (KF) reaction. This shortcoming becomes even greater as the number of samples increases.

We tested the water content in different DMSO sample quantities according to the following procedure. 30 mL of Hydranal-

Methanol Rapid, Hydranal-Methanol dry or Hydranal-CompoSolver E were introduced into the titration vessel and titrated to dryness with Hydranal-Composite 5. The sample (or sample plus water) was then added and titrated with Hydranal-Composite 5.

Results are compared in the tables below.

**Table 1. Water determination in various DMSO sample quantities**

DMSO SAMPLE VOLUME	HYDRANAL-COMPOSITE 5 CONSUMPTION	WATER QUANTITY DETECTED	WATER CONTENT DETECTED
1.0 mL	0.08 mL	0.44 mg	0.044%
2.0 mL	0.14 mL	0.77 mg	0.038%
5.0 mL	0.19 mL	1.05 mg	0.021%
10.0 mL	0.35 mL	1.93 mg	0.019%
20.0 mL	0.62 mL	3.41 mg	0.017%

**Honeywell**

HONEYWELL RESEARCH CHEMICALS PORTFOLIO  
Riedel-de Haën™ Burdick & Jackson™ Fluka™



**Table 2. Different sample quantities of DMSO added to a defined water quantity (20.0 mg)**

SAMPLE	HYDRANAL-COMPOSITE 5 CONSUMPTION	WATER QUANTITY DETECTED	RECOVERY RATE
20 mg H <sub>2</sub> O	3.81 mL	20.00 mg	100.0%
20 mg H <sub>2</sub> O + 1.0 mL DMSO	3.80 mL	19.95 mg	99.75%
20 mg H <sub>2</sub> O + 2.0 mL DMSO	3.77 mL	19.79 mg	98.95%
20 mg H <sub>2</sub> O + 5.0 mL DMSO	3.71 mL	19.48 mg	97.40%
20 mg H <sub>2</sub> O + 10.0 mL DMSO	3.49 mL	18.32 mg	91.60%
20 mg H <sub>2</sub> O + 20.0 mL DMSO	3.38 mL	17.74 mg	88.70%

Coulometric tests gave comparable findings:

- Water content detected in 1 mL DMSO: 311 ppm
- Water content detected in 20 mL DMSO: 266 ppm

Control conditions with Hydranal-Water Standard 1.0 (added water quantity: 1000 ppm) in the presence of DMSO produced the following results:

- Water content in the presence of 1 mL DMSO: 994 ppm
- Water content in the presence of 20 mL DMSO: 910 ppm

The volumetric and coulometric water content determination tests were carried out with two different samples.

The problem was picked up again in a series of subsequent tests. We ascertained that it is also impossible to perform indirect determination in a KF oven. A temperature ramp from 50°C to 250°C showed that the water is released by azeotropic distillation at temperatures between 130°C and 190°C. It is impossible to separate the water from the DMSO. The titration vessel is influenced in the same way as by direct injection of the sample. The error is comparable to the effect described for direct injection.

## Conclusion

As a result of the various facts, the recommendation must be made that no more than 1 mL of DMSO should be analyzed. The injection of several samples in a coulometric cell should be avoided. Addition of a 1000 ppm standard shows the current effect on the titration vessel.

34805 HYDRANAL-Composite 5  
34741 HYDRANAL-Methanol dry

34734 HYDRANAL-CompoSolver E  
37817 HYDRANAL-Methanol Rapid



**Europe and International**

**Thomas Wendt**

HYDRANAL Center of Excellence  
Tel: +49-5137 999-353  
Fax: +49-5137 999-698  
hydranal@honeywell.com

WATER STANDARDS

---

34828 HYDRANAL-Water Standard 1.0

34426 HYDRANAL-CRM Water Standard 1.0

AUXILIARIES

---

34241 HYDRANAL-Molecular Sieve 0.3 nm

34788 HYDRANAL-Humidity Absorber



**Europe and International**

**Agnieszka Kossakowska**

HYDRANAL Technical Specialist  
Tel: +48 512 355 628  
hydranal@honeywell.com



**USA and Canada**

**Doug Clark**

HYDRANAL Technical Center  
Tel: 1-800-Hydranal  
(1-800-493-7262)  
hydranal@honeywell.com

**To order, please contact:**

**Greyhound Chromatography and Allied Chemicals**

6 Kelvin Park  
Birkenhead, Merseyside, CH41 1LT  
Tel: 0151 649 4000  
Email: [info@greyhoundchrom.com](mailto:info@greyhoundchrom.com)  
[www.greyhoundchrom.com](http://www.greyhoundchrom.com)

**Honeywell Specialty Chemicals Seelze GmbH**

Wunstorferstrasse 40  
30926 Seelze, Germany  
Tel.: +49 (0)5137-999-353  
Fax: +49 (0)5137-999-698  
[lab-honeywell.com](http://lab-honeywell.com)

All statements and information provided herein are believed to be accurate and reliable, but are presented without guarantee, warranty or responsibility of any kind, express or implied. Statements or suggestions concerning possible use of our products are made without representation or warranty that any such use is free of patent infringement, and are not recommendations to infringe any patent. The user should not assume that all safety measures are indicated herein, or that other measures may not be required. User assumes all liability for use of the information and results obtained. WITHOUT LIMITING THE FOREGOING, HONEYWELL DISCLAIMS THE WARRANTY OF MERCHANTABILITY, FITNESS FOR USE AND NON-INFRINGEMENT. © 2018 Honeywell International Inc.



**RESPONSIBLE CARE**<sup>®</sup>  
OUR COMMITMENT TO SUSTAINABILITY

**Honeywell**

---

